

AD-A221 421

USAARL REPORT NO. 81-7

C



OPERATOR'S MANUAL FOR VARIABLE WEIGHT,
VARIABLE C.G. HELMET SIMULATOR
(Reprint)

DTIC
ELECTE
MAY 02 1990
S D OS D

DTIC FILE COPY

By
Craig M. Svoboda
James C. Warrick
Simula, Inc.

BIODYNAMICS RESEARCH DIVISION

DISTRIBUTION STATEMENT A

Approved for public release;
Distribution Unlimited

September 1981

U.S. ARMY AEROMEDICAL RESEARCH LABORATORY
FORT RUCKER, ALABAMA 36362

90 05 01 027

USAARL

NOTICE

Qualified Requesters

Qualified requesters may obtain copies from the Defense Technical Information Center, Cameron Station, Alexandria, Virginia. Orders will be expedited if placed through the librarian or other person designated to request documents from the Defense Technical Information Center.

Change of Address

Organizations receiving reports from the US Army Aeromedical Research Laboratory on automatic mailing lists should confirm correct address when corresponding about laboratory reports.

Disposition

Destroy this report when it is no longer needed. Do not return it to the originator.


Disclaimer

The views, opinions, and/or findings contained in this report are those of the authors and should not be construed as an official Department of the Army position, policy, or decision, unless so designated by other official documentation. Citation of trade names in this report does not constitute an official Department of the Army endorsement or approval of the use of such commercial items.

Reviewed:




KENT A. KIMBALL
Dir, Extramural Research Program

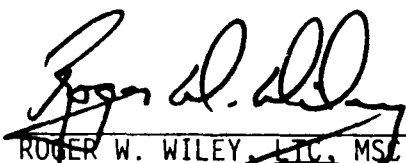

ROGER W. WILEY, LTC, MSJ
Chairman, Scientific Review
Committee

Released for Publication:

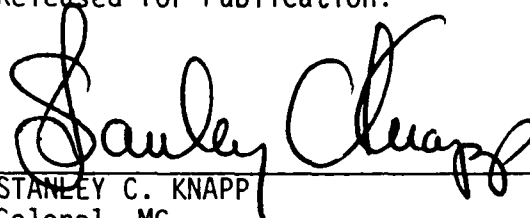

STANLEY C. KNAPP
Colonel, MC
Commanding

REPORT DOCUMENTATION PAGE		READ INSTRUCTIONS BEFORE COMPLETING FORM
1. REPORT NUMBER USAARL Rpt No. 81-7	2. GOVT ACCESSION NO.	3. RECIPIENT'S CATALOG NUMBER
4. TITLE (and Subtitle) Operator's Manual for Variable Weight, Variable C.G. Helmet Simulator		5. TYPE OF REPORT & PERIOD COVERED Reprint
		6. PERFORMING ORG. REPORT NUMBER
7. AUTHOR(s) Craig Svoboda James C. Warrick		8. CONTRACT OR GRANT NUMBER(s) DAMD-17-80-C-0089
9. PERFORMING ORGANIZATION NAME AND ADDRESS Simula Inc. 2223 South 48th Street Tempe, Arizona 85282		10. PROGRAM ELEMENT, PROJECT, TASK AREA & WORK UNIT NUMBERS
11. CONTROLLING OFFICE NAME AND ADDRESS SGRD-UAD-IV US Army Aeromedical Research Laboratory Ft. Rucker, AL 36362		12. REPORT DATE September 1981
		13. NUMBER OF PAGES 21
14. MONITORING AGENCY NAME & ADDRESS (if different from Controlling Office) USAMRDC Fort Deterick Frederick, Md 21701		15. SECURITY CLASS. (of this report) Unclassified
		15a. DECLASSIFICATION/DOWNGRADING SCHEDULE
16. DISTRIBUTION STATEMENT (of this Report) Approved for public release; distribution unlimited.		
17. DISTRIBUTION STATEMENT (of the abstract entered in Block 20, if different from Report)		
18. SUPPLEMENTARY NOTES Reprinted with permission, Simula, Inc., Tempe, Arizona, Letter dated 4 June 81		
19. KEY WORDS (Continue on reverse side if necessary and identify by block number) helmet simulator variable helmet weight variable helmet center of gravity SPH-4 Helmet <i>Continued on reverse side</i>		
20. ABSTRACT (Continue on reverse side if necessary and identify by block number) A variable weight, variable CG helmet simulator has been designed to measure the effect of US Army headgear on muscle loading and fatigue. The helmet simulator consists of two weight concealment boxes attached to opposite sides of a supporting head which in turn is mounted on the wearer's head by a suspension system taken from an SPH-4 helmet. The weight and CG can be altered by positioning variable weights within the concealment boxes. < —		

FORWARD

The helmet simulator was developed under a subcontract of U. S. Army Contract DAMD-17-80-C-0089 awarded to Wright State University, Dayton, Ohio. The subcontractor was Simula, Inc., Tempe, Arizona.

Mr. Craig M. Svoboda, project engineer, and Mr. James C. Warrick, project manager, Simula, Inc., did the developmental work and prepared this report. The U. S. Army Aeromedical Research Laboratory (USAARL) requested and received permission from Simula, Inc., to reprint and distribute the report.

CPT John C. Johnson,* Biodynamics Research Division, USAARL, served as the contracting officer's technical representative for the project.

Copies of this report have been requested by other government agencies; therefore, it is being reprinted to insure wider distribution. The contract efforts support the U. S. Army Aeromedical Research Laboratory's ongoing biodynamic research projects.



Accession For	
NTIS CRA&I	☒
DTIC TAB	☐
Unannounced	☐
Justification	
By _____	
Distribution /	
Availability Codes	
Dist	Avail and/or Special
A-1	

*CPT John C. Johnson is presently assigned to the Letterman Army Institute of Research, Persidio of San Francisco, California.

**OPERATOR'S MANUAL
FOR
VARIABLE WEIGHT,
VARIABLE C.G.
HELMET SIMULATOR**

Prepared for:

**Wright State University
Dayton, Ohio 45435**

Prepared by:

**Simula Inc.
2223 South 48th Street
Tempe, Arizona 85282**

Prepared by:

Craig M. Svoboda
**Craig M. Svoboda
Project Engineer**

Approved by:

James C. Warrick
**James C. Warrick
Project Manager**

OPERATOR'S MANUAL
FOR
VARIABLE WEIGHT, VARIABLE C.G.
HELMET SIMULATOR

1.0 DESCRIPTION

The helmet simulator consists of two weight concealment boxes attached to opposite sides of a support ring (headring) which in turn is supported upon the wearer's head by a suspension system taken from an SPH-4 helmet. The weight and c.g. can be altered by positioning variable weights within the concealment boxes. Fabric covers over the boxes prevent the test subjects from obtaining visual clues as to the c.g. location.

The minimum weight of the helmet simulator, without any variable weights in the boxes, is 2.5 lb, slightly less than the weight of most quality crash helmets made by reputable manufacturers. The addition of variable weights to the boxes can alter the center of gravity to simulate the effect of equipment attached to the outside of a helmet. The helmet simulator has been calibrated for weights of 3.2, 4.0, 5.0, 7.0, and 9.0 lb for each of the c.g. locations shown in Table 1. Figure 1 illustrates the range of c.g. variations together with definition of the coordinate axes by which the c.g. locations are measured.

As shown in Figure 1, a point midway between the left and right ear canals has been chosen as the origin of the coordinate axes. The helmet simulator has been provided with adjustment to ensure that an index point on it can be aligned with the ear canals, and also with independent adjustment to permit the suspension system to be made comfortable.

TABLE 1. SPECIFIC C.G. LOCATIONS FOR WHICH
THE HELMET SIMULATOR IS CALIBRATED

C.G. Location Description*	Displacement (cm)**		
	X	Y	Z
<u>Central</u>			
Maximum Height	0	0	8.0
Medium Height	0	0	4.0
Low Height	0	0	0
<u>Forward</u>			
Maximum Height	5.0	0	8.0
Medium Height	5.0	0	4.0
Low Height	5.0	0	0
<u>Aftward</u>			
Maximum Height	-2.5	0	8.0
Medium Height	-2.5	0	4.0
Low Height	-2.5	0	0
<u>Central</u>			
Left, Maximum Height	0	2.5	8.0
Left, Medium Height	0	2.5	4.0
Left, Low Height	0	2.5	0
<u>Central</u>			
Right, Maximum Height	0	-2.5	8.0
Right, Medium Height	0	-2.5	4.0
Right, Low Height	0	-2.5	0
<u>Forward</u>			
Left, Maximum Height	4.3	1.8	8.0
Left, Medium Height	4.3	1.8	4.0
Left, Low Height	4.3	1.8	0
<u>Forward</u>			
Right, Maximum Height	4.3	-1.8	8.0
Right, Medium Height	4.3	-1.8	4.0
Right, Low Height	4.3	-1.8	0

*C.G. locations for total weights of less than 4 lb may differ from values shown in this table. See Tables 2 through 8 for exact locations.

**Displacement from head and neck c.g. (axis directions as defined in Figure 1).

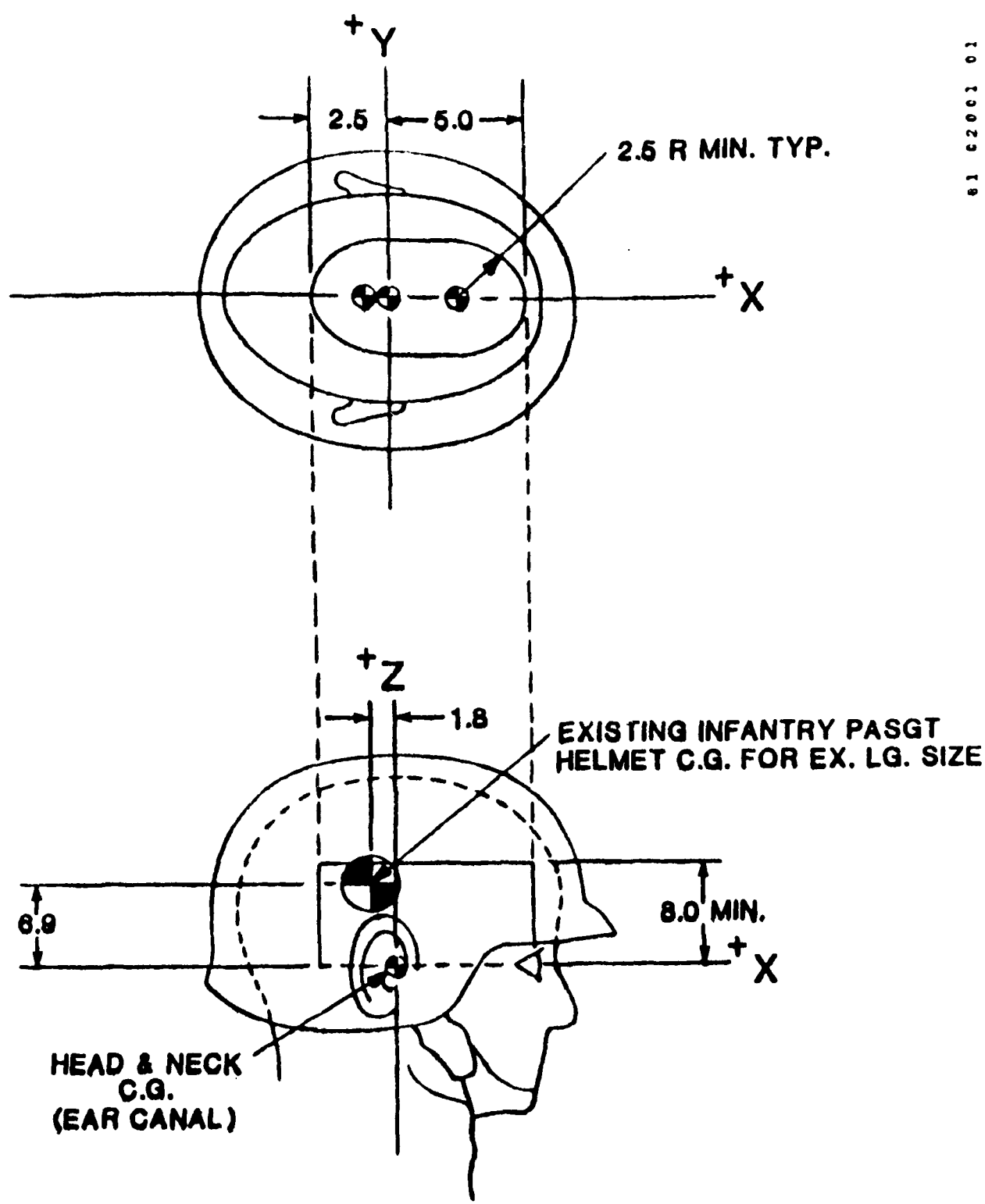


Figure 1. C.G. variation required for test helmet(s)
(all dimensions in centimeters).

2.0 ADJUSTMENT TO FIT WEARER'S HEAD

Fitting of the helmet simulator to the subject's head should be as follows:

- 2.1 Remove all weights, weight support tubes, and fabric covers from the concealment boxes.
- 2.2 Place the device on the subject's head.
- 2.3 Adjust the headband suspension buckle.
- 2.4 While applying downward pressure upon the helmet handle to compress the hair, adjust the crown strap buckles (3 places) to permit comfortable vertical positioning of the earcups on the wearer's ears.
- 2.5 Snap and adjust chin strap, also adjust rear earcup retainer strap buckle.
- 2.6 With helmet sitting squarely upon subject's head, readjust all straps.
- 2.7 Determine if ear canals are aligned with holes in concealment boxes. If not, loosen the screw attaching one of the weight boxes to the head ring, slide the box to the correct position, then retighten the screw. Repeat for the other box.
- 2.8 Repeat the preceding steps, if necessary, to obtain snug and comfortable fit.

3.0 ADJUSTMENT OF VARIABLE WEIGHTS

Weights can be installed in the concealment boxes, either with the helmet in place on the subject's head, or with the helmet removed and placed upon a table out of view of the subject.


Tables 2 through 8 identify, for chosen combinations of weights and c.g. locations, the quantity of variable weights which must be added to each concealment box and the location of the weights within each box.

The installation of weights can best be understood by studying the following example:

EXAMPLE: Suppose that it has been chosen to test a c.g. location that is forward, to the left, and high. This condition is found in Table 7. Now suppose the total desired weight is 4.0 lb. This condition is found in row two of that table.

Remove the weight tube from the left-hand weight box (if this has not already been done) by pressing on the button and swinging the top of the tube out of the box. Pull the lock pin in the master weight out of engagement with the tube but not completely out of the master weight. Reposition the master weight on the tube so that graduation 22 appears in the window formed by the notch in the master weight as shown in Figure 2, and re-engage the lock pin with the tube. Additional weights, one each 0.2 lb and 0.5 lb, should then be assembled to that side of the master weight which has higher numbered tube graduations.

After loading the left-hand tube with the correct weights, install it by inserting the 3/16-in. diameter end of the tube into hole V in the bottom of the left-hand box. Depress the button at the top of the tube and swing the top of the tube into the box, allowing the button to engage with that hole which is in line with the bottom of the tube.

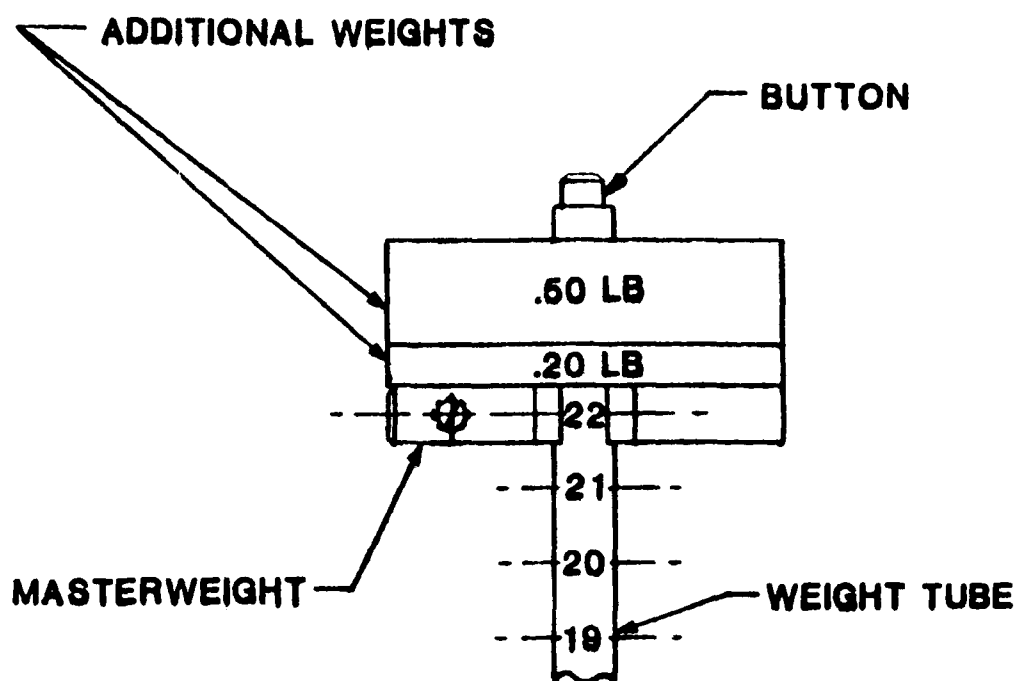

61 02002 02

Figure 2. Appearance of weights loaded onto weight tube.

Verify correctness of weight location and quantity. Notice that both the left-hand weight tube and weight concealment box are color coded yellow.

Attach fabric cover, using velcro patches.

Follow similar procedure for the right-hand box: position the master weight at graduation 22; add one each weights of 0.1 and 0.2 lb; and position tube in right-hand box at hole U. Verify correctness, noting that both right-hand weight tube and right-hand box are color coded white. It is NOT necessary for left-hand and right-hand weights to be positioned at identical graduations.

4.0 RECALIBRATION OF HELMET SIMULATOR

Refer to Appendix A for information concerning recalibration if it becomes necessary to perform tests with weight and c.g. combinations other than those for which the helmet simulator has already been calibrated.

TABLE 2. CENTRAL X, CENTRAL Y, VARIABLE Z, AND WEIGHT

C.G. Location (cm)			Total Weight (lb)	(Yellow color code) Left-hand (L.H.) Box							Right-hand (R.H.) Box						
				Position of Master Weight on L.H. Tube	Additional Weights (lb)				Position of L.H. Tube in L.H. Box	Position of Master Weight on R.H. Tube	Additional Weights (lb)				Position of R.H. Tube in R.H. Box		
					0.1	0.2	0.5	1.0			0.1	0.2	0.5	1.0			
																Quantity	Quantity
0	0	6.8	3.2	23	H	1	-	-	-	-	23	1	-	-	-	H	
0	0	8	4	22	J	-	-	1	-	-	22	-	-	1	-	J	
		High	5	19	J	-	-	-	1	-	19	-	-	-	1	J	
			7	16	K	-	-	-	2	-	16	-	-	-	2	K	
			9	14	K	-	-	-	3	-	14	-	-	-	3	K	
4			3.2	9	H	1	-	-	-	-	9	1	-	-	-	H	
		Med	4	11	J	-	-	1	-	-	11	-	-	1	-	J	
			5	10	J	-	-	-	1	-	10	-	-	-	1	J	
			7	10	K	-	-	-	2	-	10	-	-	-	2	K	
			9	8	K	-	-	-	3	-	8	-	-	-	3	K	
2.6			3.2	2	H	1	-	-	-	-	2	1	-	-	-	H	
1			4	2	J	-	-	1	-	-	2	-	-	1	-	J	
0			5	2	J	-	-	-	1	-	2	-	-	-	1	J	
0			7	3	K	-	-	-	2	-	3	-	-	-	2	K	
0			9	3	K	-	-	-	3	-	3	-	-	-	3	K	

TABLE 3. FORWARD X, CENTRAL Y, VARIABLE Z, AND WEIGHT

C.G. Location (cm)			Total Weight (lb)	(Yellow color code) Left-hand (L.H.) Box						Right-hand (R.H.) Box									
				Position of Master Weight on L.H. Tube	Additional Weights (lb)				Position of L.H. Tube in L.H. Box	Position of Master Weight on R.H. Tube	Additional Weights (lb)				Position of R.H. Tube in R.H. Box				
					0.1	0.2	0.5	1.0			0.1	0.2	0.5	1.0					
																Quantity			
X	Y	Z	4	0	6.8	3.2	23	1	-	-	-	-	23	1	-	-	-	-	Z
5	0	8	4			22	22	-	-	1	-	-	22	-	-	1	-	-	X
		High	5			19	19	-	-	-	1	-	19	-	-	-	1	-	T
			7			16	16	-	-	-	2	-	16	-	-	-	2	-	R
			9			14	14	-	-	-	3	-	14	-	-	-	3	-	R
4	0	4	3.2			9	9	1	-	-	-	-	9	1	-	-	-	-	Z
5	0	4	4			11	11	-	-	1	-	-	11	-	-	1	-	-	X
		Med	5			10	10	-	-	-	1	-	10	-	-	-	1	-	T
			7			10	10	-	-	-	2	-	10	-	-	-	2	-	R
			9			8	8	-	-	-	3	-	8	-	-	-	3	-	R
4	0	2.6	3.2			2	2	1	-	-	-	-	2	1	-	-	-	-	Z
5	0	1	4			2	2	-	-	1	-	-	2	-	-	1	-	-	X
		0	5			2	2	-	-	-	1	-	2	-	-	-	1	-	T
		0	7			3	3	-	-	-	2	-	3	-	-	-	2	-	R
		0	9			3	3	-	-	-	3	-	3	-	-	-	3	-	R

TABLE 4. AFTWARD X, CENTRAL Y, VARIABLE Z, AND WEIGHT

C.G. Location (cm)			Total Weight (lb)	(Yellow color code)						Right-hand (R.H.) Box											
				Left-hand (L.H.) Box						Position of Master Weight on R.H. Tube	Additional Weights (lb)			Position of Master Weight on R.H. Tube	Additional Weights (lb)			Position of R.H. Tube in R.H. Box			
				Position of Master Weight on L.H. Tube	Additional Weights (lb)			Position of L.H. Tube in L.H. Box	Quantity		0.1	0.2	0.5		1.0	Quantity	0.1		0.2	0.5	1.0
					0.1	0.2	0.5														
-1.4	0	6.8	3.2	23	1	-	-	-	A	23	1	-	-	-	-	-	-	A			
-2.5	0	8	4	22	-	-	1	-	C	22	-	-	1	-	-	-	-	C			
		High	5	19	-	-	-	1	E	19	-	-	-	-	1	-	-	E			
			7	16	-	-	-	2	F	16	-	-	-	-	2	-	-	F			
			9	14	-	-	-	3	G	14	-	-	-	-	3	-	-	G			
-1.4	0	4	3.2	9	1	-	-	-	A	9	1	-	-	-	-	-	-	A			
-2.5	0	4	4	11	-	-	1	-	C	11	-	-	1	-	-	-	-	C			
		Med	5	11	-	-	-	1	E	11	-	-	-	-	1	-	-	E			
			7	10	-	-	-	2	F	10	-	-	-	-	2	-	-	F			
			9	8	-	-	-	3	G	8	-	-	-	-	3	-	-	G			
-1.4	0	2.6	3.2	2	1	-	-	-	A	2	1	-	-	-	-	-	-	A			
-2.5	0	1	4	2	-	-	1	-	C	2	-	-	1	-	-	-	-	C			
		0	5	2	-	-	-	1	E	2	-	-	-	-	1	-	-	E			
		0	7	3	-	-	-	2	F	3	-	-	-	-	2	-	-	F			
		0	9	3	-	-	-	3	G	3	-	-	-	-	3	-	-	G			

TABLE 5. CENTRAL X, LEFT Y, VARIABLE Z, AND WEIGHT

C.G. Location (cm)			Total Weight (lb)	(Yellow color code) Left-hand (L.H.) Box						Right-hand (R.H.) Box					
				Position of Master Weight on L.H. Tube	Additional Weights (lb)				Position of L.H. Tube in L.H. Box	Position of Master Weight on R.H. Tube	Additional Weights (lb)				Position of R.H. Tube in R.H. Box
					0.1	0.2	0.5	1.0			0.1	0.2	0.5	1.0	
X	Y	Z													
0	1.3	6.8	3.2	23	-	1	-	-	H	23	-	-	-	H	
0	2.5	8	4	22	1	1	1	-	J	21	-	1	-	J	
			5	19	1	1	-	1	J	18	-	1	-	J	
		Hgh	7	16	-	-	1	2	J	16	-	-	1	J	
			9	14	1	-	1	3	J	14	-	2	-	J	
0	1.3	4	3.2	9	2	-	-	-	H	9	-	-	-	F	
0	2.5	4	4	11	1	1	1	-	J	10	-	1	-	J	
		Med	5	11	1	1	-	1	J	10	-	1	-	J	
			7	9	-	-	1	2	J	9	-	-	1	J	
			9	8	1	-	1	3	J	8	-	2	-	J	
0	1.3	2.6	3.2	2	-	1	-	-	H	2	-	-	-	H	
0	2.5	1	4	2	1	1	1	-	J	2	-	1	-	J	
		0	5	2	1	1	-	1	J	2	-	1	-	J	
		0	7	3	-	-	1	2	J	3	-	-	1	J	
		0	9	3	1	-	1	3	J	3	-	2	-	J	

TABLE 6. CENTRAL X, RIGHT Y, VARIABLE Z, AND WEIGHT

C.G. Location (cm.)			Total Weight (lb)	(Yellow color code) Left-hand (L.H.) Box							Right-hand (R.H.) Box						
				Position of Master Weight on L.H. Tube	Additional Weights (lb)				Position of L.H. Tube in L.H. Box	Position of Master Weight on R.H. Tube	Additional Weights (lb)				Position of R.H. Tube in R.H. Box		
					Quantity						Quantity						
					0.1	0.2	0.5	1.0			0.1	0.2	0.5	1.0			
0	-1.3	6.8	3.2	23	-	-	-	-	H	23	-	1	-	-	H		
0	-2.5	8	4	21	-	1	-	-	J	22	1	1	1	-	J		
			5	18	-	1	1	-	J	19	1	1	-	1	J		
		High	7	16	-	-	1	1	J	16	-	-	1	2	J		
			9	14	-	2	-	2	J	14	1	-	1	3	J		
0	-1.3	4	3.2	9	-	-	-	-	H	9	-	1	-	-	H		
0	-2.5	4	4	10	-	1	-	-	J	11	1	1	1	-	J		
		Med	5	10	-	1	1	-	J	11	1	1	-	1	J		
			7	9	-	-	1	1	J	9	-	-	1	2	J		
			9	8	-	2	-	2	J	8	1	-	1	3	J		
0	-1.3	2.6	3.2	2	-	-	-	-	H	2	-	1	-	-	H		
0	-2.5	1	4	2	-	1	-	-	J	2	1	1	1	-	J		
		0	5	2	-	1	1	-	J	2	1	1	-	1	J		
		0	7	3	-	-	1	1	J	3	-	-	1	2	J		
		0	9	3	-	2	-	2	J	3	1	-	1	3	J		

TABLE 7. FORWARD X, LEFT Y, VARIABLE Z, AND WEIGHT

C.G. Location (cm)			Total Weight (lb)	(Yellow color code)							Right-hand (R.H.) Box								
				Left-hand (L.H.) Box							Position of Master Weight on R.H. Tube	Additional Weights (lb)			Position of R.H. Tube in F.H. Box				
				Position of Master Weight on L.H. Tube	Additional Weights (lb)			Position of L.H. Tube in L.H. Box	Quantity										
X	Y	Z								0.1	0.2	0.5	1.0		0.1	0.2	0.5	1.0	
4.0	1.3	6.8	3.2	23	-	1	-	-	Z	-	-	-	-	-	-	-	-	-	Z
4.3	1.8	8	4	22	-	1	1	-	V	-	1	-	-	1	-	-	-	-	U
		High	5	19	-	1	-	1	S	-	-	1	-	1	1	-	-	-	S
			7	16	1	1	-	2	R	-	-	-	1	1	1	1	1	1	Q
			9	14	-	2	-	3	Q	-	-	-	-	-	-	1	2	2	P
4.0	1.3	4	3.2	9	-	1	-	-	Z	-	-	-	-	-	-	-	-	-	Z
4.3	1.8		4	10	-	1	1	-	V	-	1	-	-	1	-	-	-	-	U
		Med	5	10	-	1	-	1	S	-	-	1	-	1	1	1	-	-	S
			7	9	1	1	-	2	R	-	-	-	1	1	1	1	1	1	Q
			9	8	-	2	-	3	Q	-	-	-	-	-	-	1	2	2	P
4.0	1.3	2.6	3.2	2	-	1	-	-	Z	-	-	-	-	-	-	-	-	-	Z
4.3	1.8	1	4	2	-	1	1	-	V	-	1	-	-	1	-	-	-	-	U
		0	5	2	-	1	-	1	S	-	-	1	-	1	1	1	-	-	S
		0	7	4	1	1	-	2	R	-	-	-	1	1	1	1	1	1	Q
		0	9	3	-	2	-	3	Q	-	-	-	-	-	-	1	2	2	P

TABLE 8. FORWARD N, RIGHT Y, VARIABLE Z, AND WEIGHT

C.G. Location (cm)			Total Weight (lb)	(Yellow color code) Left-hand (L.H.) Box							Right-hand (R.H.) Box								
				Position of Master Weight on L.H. Tube	Additional Weights (lb)				Position of L.H. Tube in L.H. Box	Position of Master Weight on R.H. Tube	Additional Weights (lb)				Position of R.H. Tube in R.H. Box				
					0.1	0.2	0.5	1.0			0.1	0.2	0.5	1.0					
																Quantity			
X	Y	Z																	
4.0	-1.3	6.8	3.2	23	-	-	-	-	Z	23	-	1	-	-	Z				
4.3	-1.8	8	4	22	1	1	-	-	U	22	-	1	1	-	V				
		High	5	19	1	1	1	-	S	19	-	1	-	1	S				
			7	16	-	1	1	1	Q	16	1	1	-	2	R				
			9	14	1	-	1	2	P	14	-	2	-	3	Q				
4.0	-1.3	4	3.2	9	-	-	-	-	Z	9	-	1	-	-	Z				
4.3	-1.8		4	11	1	1	-	-	U	10	-	1	1	-	V				
		Med	5	10	1	1	1	-	S	10	-	1	-	1	S				
			7	10	-	1	1	1	Q	9	1	1	-	2	R				
			9	8	1	-	1	2	P	8	-	2	-	3	Q				
4.0	-1.3	2.6	3.2	2	-	-	-	-	Z	2	-	1	-	-					
4.3	-1.8	1	4	2	1	1	-	-	U	2	-	1	1	-	V				
		0	5	2	1	1	1	-	S	2	-	1	-	1	S				
		0	7	3	-	1	1	1	Q	4	1	1	-	2	R				
		0	9	3	1	-	1	2	P	3	-	2	-	3	Q				

APPENDIX A

RECALIBRATION

1. Place helmet on scale and add variable weights to provide the desired total weight.
2. Assemble the balsa wood calibration brackets to the helmet and drill into each bracket a hole which is in line with the desired c.g. (see Figure A-1).
3. Hang the helmet from a string or cable passing through the new hole in one bracket, and shift the weights to level the helmet. (A bubble level attached to the helmet may facilitate this.) See Figure A-2.
4. Repeat for the hole in the other bracket (See Figure A-3).
5. Record the weight quantity and location on a copy of data sheet (Figure A-4).

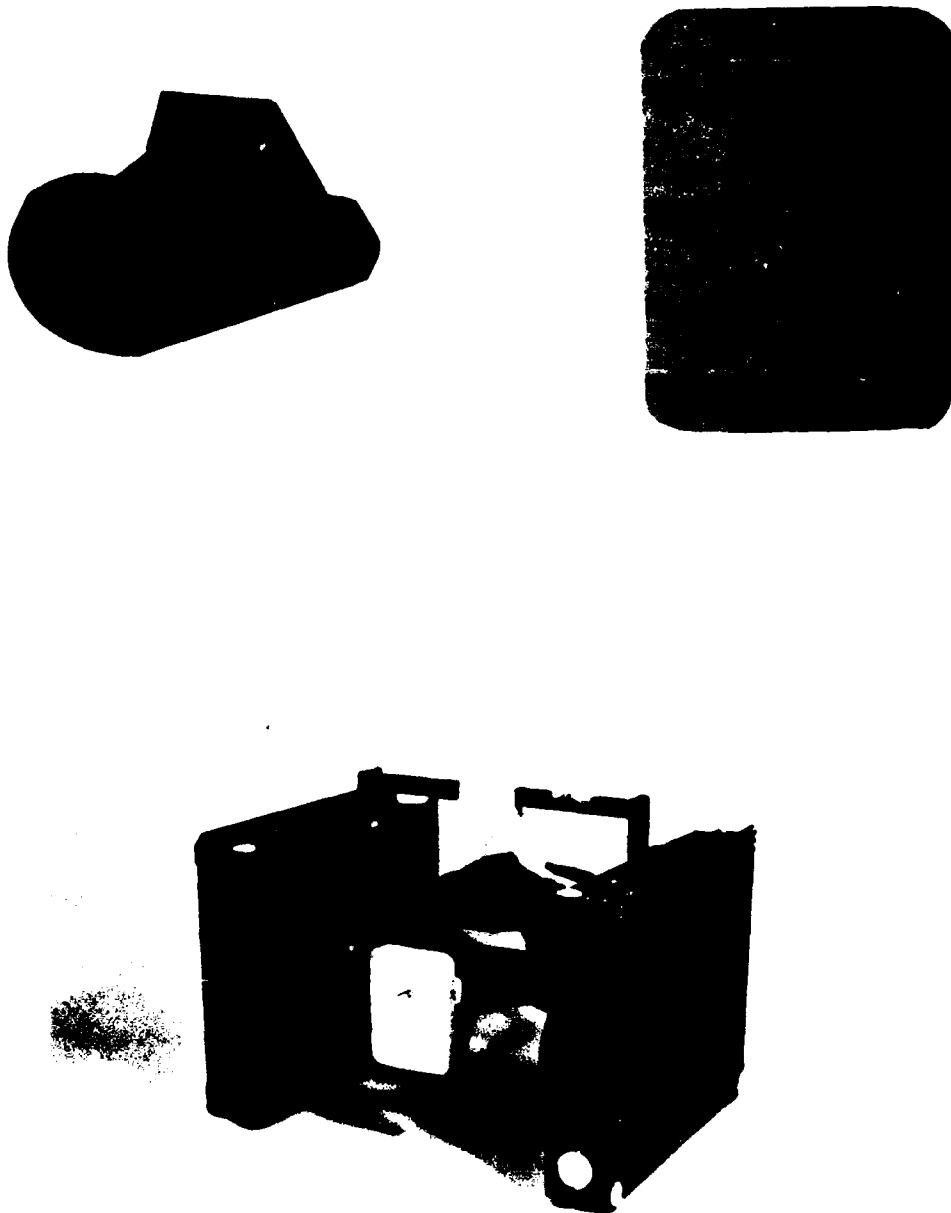


Figure A-1. Balsa wood calibration brackets and assembly on helmet.



Figure A-2. Suspend helmet by cable attached to proper hole in top bracket. Shift weights between left-hand and right-hand boxes to balance laterally, and shift weight tubes from hole-to-hole to balance in the fore-and-aft direction.


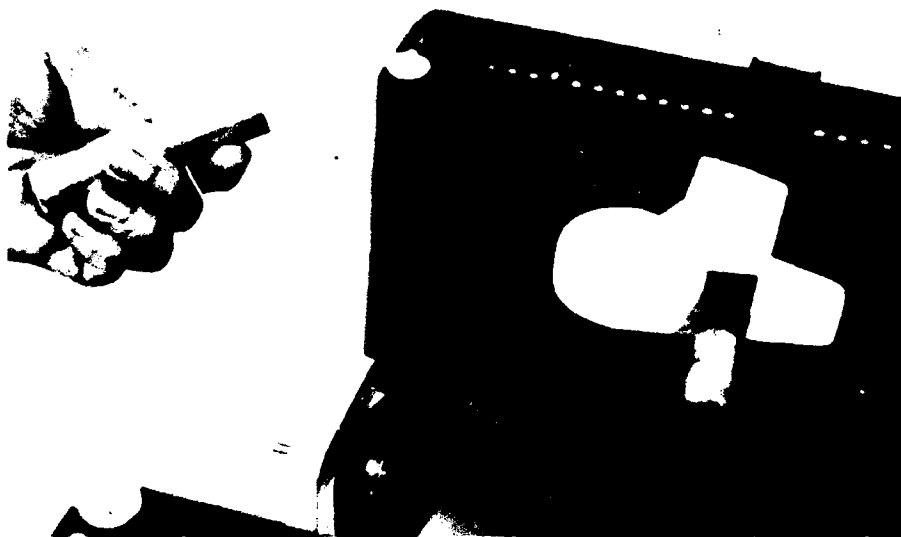


Figure A-3. Relocate cable to proper hole in front bracket, hang helmet and adjust position of weights on each weight tube to balance helmet with aid of bubble level.

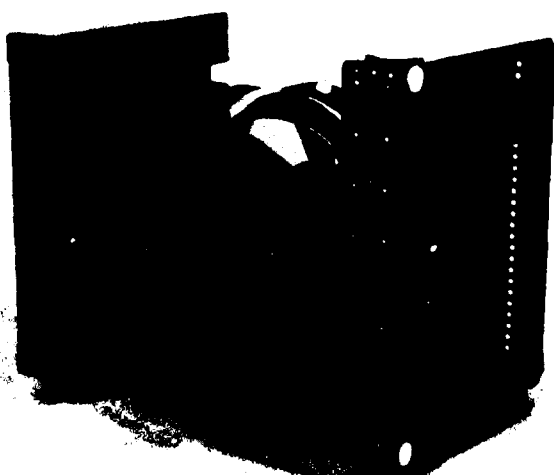

C.G. Location (cm.)			Total Weight (lb)	Left-hand (L.H.) Box					Right-hand (R.H.) Box									
				Position of Master Weight on L.H. Tube	Additional Weights (lb)				Position of L.H. Tube in L.H. Box	Position of Master Weight on R.H. Tube	Additional Weights (lb)				Position of R.H. Tube in R.H. Box			
					0.1	0.2	0.5	1.0			0.1	0.2	0.5	1.0				
X	Y	Z						Quantity					Quantity					
		High	4															
			5															
			7															
			9															
		Med	4															
			5															
			7															
			9															
		Low	4															
			5															
			7															
			9															

Figure A-4. Sample data sheet.

INITIAL DISTRIBUTION

Defense Technical Information Center Cameron Station Alexandria, VA 22314 (12)	Aeromechanics Laboratory US Army Research & Technology Labs Ames Research Center, M/S 215-1 Moffett Field, CA 94035 (1)
Under Secretary of Defense for Research and Engineering ATTN: Military Assistant for Medical and Life Sciences Washington, DC 20301 (1)	Sixth United States Army ATTN: SMA Presidio of San Francisco, California 94129 (1)
Uniformed Services University of the Health Sciences 4301 Jones Bridge Road Bethesda, MD 20014 (1)	Director Army Audiology & Speech Center Walter Reed Army Medical Center Forest Glen Section, Bldg 156 Washington, DC 20012 (1)
Commander US Army Medical Research and Development Command ATTN: SGRD-RMS/Ms. Madigan Fort Detrick Frederick, MD 21701 (5)	Harry Diamond Laboratories Scientific & Technical Information Offices 2800 Powder Mill Road Adelphi, MD 20783 (1)
Redstone Scientific Information Center ATTN: DRDMI-TBD US Army Missile R&D Command Redstone Arsenal, AL 35809 (1)	US Army Ordnance Center & School Library, Bldg 3071 ATTN: ATSL-DOSL Aberdeen Proving Ground, MD 21005 (1)
US Army Yuma Proving Ground Technical Library Yuma, AZ 85364 (1)	US Army Environmental Hygiene Agency Library, Bldg E2100 Aberdeen Proving Ground, MD 21010 (1)
US Army Aviation Engineering Flight Activity ATTN: DAVTE-M (Technical Library) Edwards AFB, CA 93523 (1)	Technical Library Chemical Systems Laboratory Aberdeen Proving Ground, MD 21010 (1)
US Army Combat Developments Experimentation Command Technical Library HQ, USACDEC Box 22 Fort Ord, CA 93941 (1)	US Army Materiel Systems Analysis Agency ATTN: Reports Distribution Aberdeen Proving Ground, MD 21005 (1)

Commander
US Army Medical Research Institute
of Chemical Defense
Aberdeen Proving Ground
21010

(1)

HQ, First United States Army
ATTN: AFKA-MD (Surgeon's Office)
Fort George G. Meade, MD 20755

(1)

Director
Ballistic Research Laboratory
ATTN: DRDAR-TSB-S (STIFF)
Aberdeen Proving Ground, MD
21005

(2)

US Army Research & Development
Technical Support Activity
Fort Monmouth, NJ 07703

(1)

Commander/Director
US Army Combat Surveillance &
Target Acquisition Laboratory
ATTN: DELCS-D
Fort Monmouth, NJ 07703

(1)

US Army Avionics R&D Activity
ATTN: DAVAA-O
Fort Monmouth, NJ 07703

(1)

US Army White Sands Missile Range
Technical Library Division
White Sands Missile Range
New Mexico 88002

(1)

Chief
Benet Weapons Laboratory
LCWSL, USA ARRADCOM
ATTN: DRDAR-LCB-TL
Watervliet Arsenal
Watervliet, NY 12189

(1)

US Army Research & Technology Labs
Propulsion Laboratory MS 100-5
NASA Lewis Research Center
Cleveland, OH 44135

(1)

US Army Field Artillery School
Library
Snow Hall, Room 16
Fort Sill, OK 73503

(1)

US Army Dugway Proving Ground
Technical Library
Bldg 5330
Dugway, UT 84022

(1)

US Army Materiel Development &
Readiness Command
ATTN: DRCSG
5001 Eisenhower Avenue
Alexandria, VA 22333

(1)

US Army Foreign Science &
Technology Center
ATTN: DRXST-IS1
220 7th St., NE
Charlottesville, VA 22901

(1)

Commander
US Army Training and Doctrine Command
ATTN: ATCD
Fort Monroe, VA 23651

(2)

Commander
US Army Training and Doctrine Command
ATTN: Surgeon
Fort Monroe, VA 23651

(1)

US Army Research & Technology Labs
Structures Laboratory Library
NASA Langley Research Center
Mail Stop 266
Hampton, VA 23665

(1)

Commander
10th Medical Laboratory
ATTN: DEHE (Audiologist)
APO New York 09180

(1)

Commander
US Army Natick R&D Laboratories
ATTN: Technical Librarian
Natick, MA 01750

(1)

Commander US Army Troop Support & Aviation Materiel Readiness Command ATTN: DRSTS-W St. Louis, MO 63102 (1)	US Air Force Armament Development & Test Center Technical Library Eglin AFB, FL 32542 (1)
Commander US Army Aviation R&D Command ATTN: DRDAV-E P. O. Box 209 St. Louis, MO 63166 (1)	US Air Force Institute of Technology (AFIT/LDE) Bldg 640, Area B Wright-Patterson AFB, OH 45433 (1)
Director US Army Human Engineering Laboratory ATTN: Technical Library Aberdeen Proving Ground, MD 21005 (1)	US Air Force Aerospace Medical Division School of Aerospace Medicine Aeromedical Library/TSK-4 Brooks AFB, TX 78235 (1)
Commander US Army Aviation R&D Command ATTN: Library P. O. Box 209 St. Louis, MO 63166 (1)	Director of Professional Services Office of The Surgeon General Department of the Air Force Washington, DC 20314 (1)
Commander US Army Health Services Command ATTN: Library Fort Sam Houston, TX 78234 (1)	Human Engineering Division Air Force Aerospace Medical Research Laboratory ATTN: Technical Librarian Wright-Patterson AFB, OH 45433 (1)
Commandant US Army Academy of Health Sciences ATTN: Library Fort Sam Houston, TX 78234 (1)	US Navy Naval Weapons Center Technical Library Division Code 2333 China Lake, CA 93555 (1)
Commander US Army Airmobility Laboratory ATTN: Library Fort Eustis, VA 23604 (1)	US Navy Naval Aerospace Medical Institute Library Bldg 1953, Code 012 Pensacola, FL 32508 (1)
Air University Library (AUL/LSE) Maxwell AFB, AL 36112 (1)	US Navy Naval Submarine Medical Research Lab Medical Library, Naval Submarine Base Box 900 Groton, CT 06340 (1)
US Air Force Flight Test Center Technical Library, Stop 238 Edwards AFB, CA 93523 (1)	

Director
Naval Biosciences Laboratory
Naval Supply Center, Bldg 844
Oakland, CA 94625

(1)

Naval Air Systems Command
Technical Library AIR 950D
Rm 278 Jefferson Plaza II
Department of the Navy
Washington, DC 20361

(1)

US Navy
Naval Research Laboratory Library
Code 1433
Washington, DC 20375

(1)

US Navy
Naval Air Development Center
Technical Information Division
Technical Support Department
Warminster, PA 18974

(1)

Human Factors Engineering Division
Aircraft & Crew Systems Technology
Directorate

Naval Air Development Center
Warminster, PA 18974

(1)

US Navy
Naval Research Laboratory Library
Shock & Vibration Information Center
Code 8404
Washington, DC 20375

(1)

Director of Biological & Medical
Sciences Division
Office of Naval Research
800 N. Quincy Street
Arlington, VA 22217

(1)

Commanding Officer
Naval Medical R&D Command
National Naval Medical Center
Bethesda, MD 20014

(1)

Commanding Officer
Naval Biodynamics Laboratory
P. O. Box 29407
New Orleans, LA 70189

(1)

FAA Civil Aeromedical Institute
ATTN: Library
Box 25082
Oklahoma City, OK 73125

(1)

Department of Defence
R.A.N. Research Laboratory
P. O. Box 706
Darlinghurst, N.S.W. 2010
Australia

(1)

Canadian Society of Avn Med
c/o Academy of Medicine, Toronto
ATTN: Ms. Carmen King
288 Bloor Street West
Toronto, Ontario
M5S 1V8

(1)

COL F. Cadigan
DAO-AMLOUS B
Box 36, US Embassy
FPO New York 09510

(1)

DCIEM/SOAM
MAJ J. Soutendam (Ret.)
1133 Sheppard Avenue West
P. O. Box 2000
Downsview, Ontario
M3M 3B9

(1)

Dr. E. Hendler
Code 6003
Naval Air Development Center
Warminster, PA 18974

(1)

Commander
US Army Transportation School
ATTN: ATSP-TD-ST
Fort Eustis, VA 23604

(1)

Staff Officer, Aerospace Medicine
RAF Staff
British Embassy
3100 Massachusetts Avenue, N.W.
Washington, DC 20008 (1)